

# RECOMMENDED CLEANROOM GOWNING PROCEDURE

Before gowning, all jewellery and make-up should be removed.  
Tacky Mats and Shoe Cleaner may also be used.

Note: These are gowning procedure examples. Each customer and facility may have a different gowning procedure. Ask questions to determine the gowning procedure at each facility.

## Suggested Nonsterile Cleanroom Gowning Procedure

### Step 1: Pre-Entry

#### A. Bouffant/Hair Net

Be sure to contain all hair.



#### B. Shoe Covers

Contain all laces and tassels.



### Step 2: Gowning

#### A. Gowning Gloves

Recommended for Class 100 and better.



#### B. Mask

Worn under hood.



#### C. Hood

Ensure snug fit and proper face/neck seal.



### Step 3: Gowning

#### Coverall

Step into the coverall. Be sure sleeves and upper garment do not touch bench or floor. Tuck in hood before zipping.



### Step 4: Gowning

#### Boot Covers

Put on boots and pull boots (hightop shoe covers) over legs of coverall.



### Step 5: Gowning

#### Goggles

Wear goggles/safety glasses when eye protection or additional particulate control is desired.



### Step 6: Gowning

#### Second set of Gloves (optional)

Place hem of glove over cuff of sleeve.



## Suggested Sterile Cleanroom Gowning Procedure

### Step 1: Pre-Entry

#### A. Bouffant/Hair Net

Be sure to contain all hair.



#### B. Shoe Covers

Contain all laces and tassels.



### Step 2: Gowning

A. Wash hands using bactericidal soap and dry thoroughly.



#### B. First Sterile Gloves

Only touch inside of glove.



#### C. Sanitise gloves after donning each article.



### Step 3: Gowning

#### A. Mask

Worn under hood.



#### B. Hood

Ensure proper neck seal.




### Step 4: Gowning

#### Coverall

Carefully remove coverall from package by holding inside neck area or seam, unfold garment while preventing it from touching you, the floor, your clothes or bench.

When inserting leg, ensure donning snaps (if used) release and foot clears elastic ankle prior to transferring.

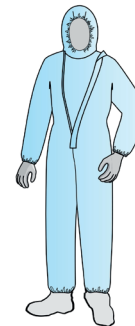

### Step 5: Gowning

#### A. Boot Covers

Put on boots and pull boots (high-top shoe covers) over legs of coverall. The first boot donned can now be placed in the "clean" side of gowning area.



B. Don second boot and place other foot in the "clean" side of the gowning area.



### Step 6: Gowning


#### A. Goggles

Adjust for a snug facial fit.




#### B. Second Sterile Gloves

Don touching only inside of gloves, placing over first pair making sure cuffs are secured over sleeves.

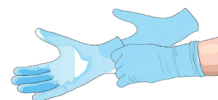

C. Perform an inspection in mirror before entering the sterile cleanroom.

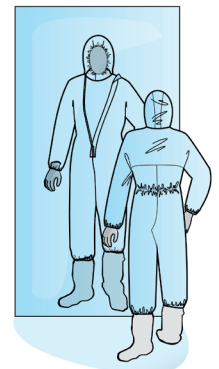

Outer Gloves



Goggles



Hood



Mask



Boot Covers



Coverall



Inner Gloves



Bouffant/Hair Net



Shoe Covers

© 2015 Thermo Fisher Scientific Inc. All rights reserved.

Trademarks used are owned as indicated at [www.fishersci.com/trademarks](http://www.fishersci.com/trademarks).

Austria: +43(0)800-20 88 40 Belgium: +32 (0)56 260 260 Denmark: +45 70 27 99 20  
Germany: +49 (0)2304 9325 Ireland: +353 (0)1 885 5854 Italy: +39 02 950 59 478  
Finland: +358 (0)9 8027 6280 France: +33 (0)3 88 67 14 14 Netherlands: +31 (0)20 487 70 00  
Norway: +47 22 95 59 59 Portugal: +351 21 425 33 50 Spain: +34 902 239 303  
Sweden: +46 31 352 32 00 Switzerland: +41 (0)56 618 41 11 UK: +44 (0)1509 555 500

 **Fisher Scientific**

A Thermo Fisher Scientific Brand